


excillum

REDEFINING THE X-RAY TUBE

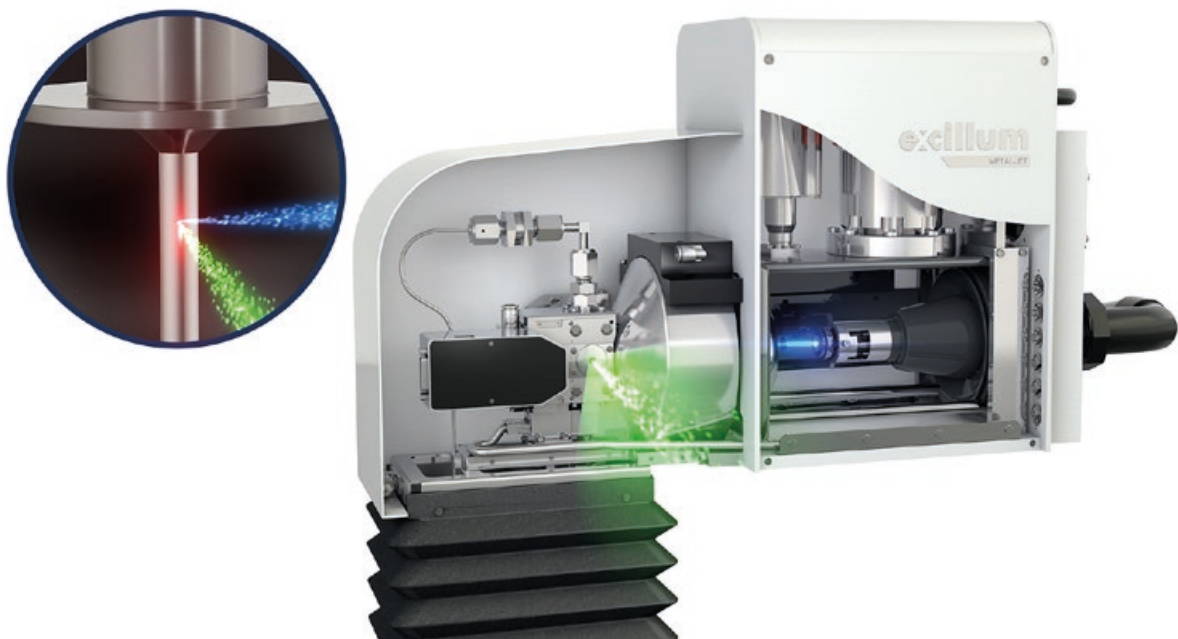

MetalJet Technology


A brief introduction

The MetalJet Concept


A conventional microfocus X-ray tube with the solid-metal anode replaced by a liquid-metal jet. The metal jet supports higher electron-beam power and can therefore generate higher X-ray flux.

Extreme performance for
5-40 μm source diameter

E-beam power density

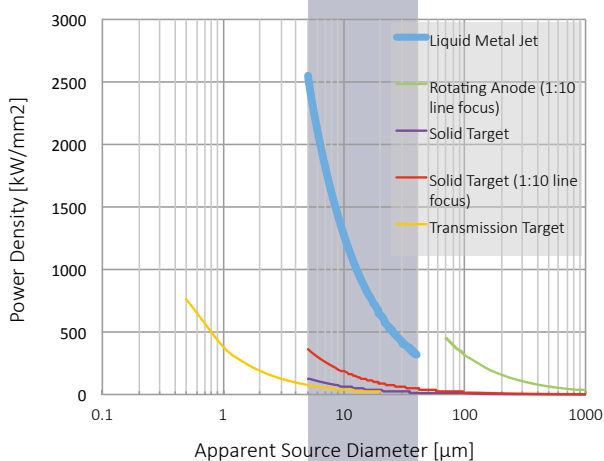

E-beam power

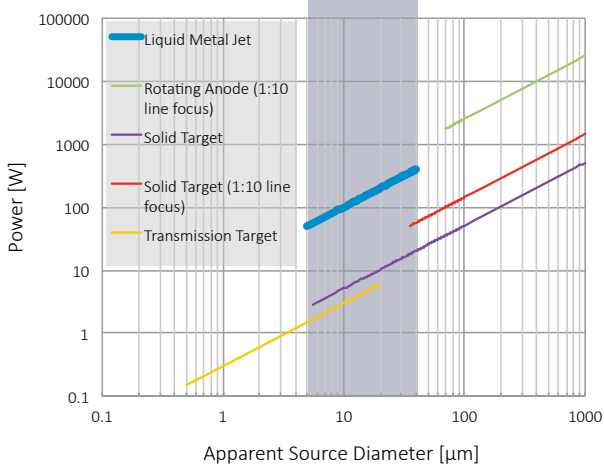

Power Loading Capability

The X-ray power of all electron-impact X-ray generators is limited by the thermal power loading of the anode. In conventional solid anode technology, the surface temperature of the anode must be well below the melting point in order to avoid damage, and this is fundamentally limited by the anode target material properties, primarily the melting point and the thermal conductivity.

The liquid-metal anode is different since the limitation to maintain the target at well below melting point is removed. This is due to the fact that the material is already molten and that it is re-generative by nature. Fresh target material is supplied at a rate of close to 100 m/s. As the anode is continuously re-generated, the electron beam and anode interaction may actually be destructive.

Extreme Brightness

Somewhat counter-intuitive, the power loading capability of small-focus X-ray tubes roughly scales with the diameter and not the area of the e-beam focus. Therefore, the brightness is inversely proportional to the source diameter.

By combining extreme power loading capability and a small electron focus, a liquid-jet source can achieve unprecedented brightness at micron spot sizes.

X-Ray Spectra of Liquid Metal

In order to reach different X-ray emission lines, different metal alloys are used. First generation metal-jet sources feature metal alloys that are molten at more or less room temperature. Still, several alloys have emission characteristics similar to regular solid anodes. Future upgrades can also include alloys with higher melting points.


Gallium Alloy

A gallium (Ga) rich alloy is available. Its 9.2 keV (1.3 Å) K α emission line is close to the copper (Cu) K α emission line.


Indium Alloy

An indium (In) rich alloy is also available. Its 24 keV (0.51 Å) K α emission line is close to the silver (Ag) K α emission line.

Spectra of Gallium alloy and Copper

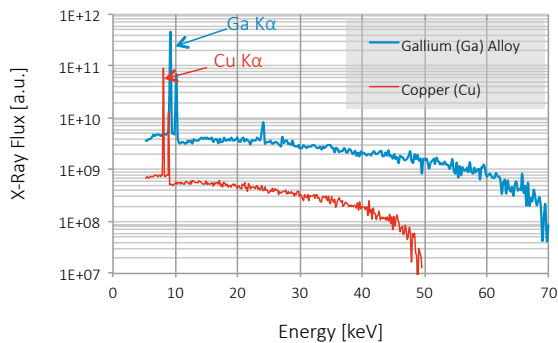


Spectra of Indium alloy and Silver

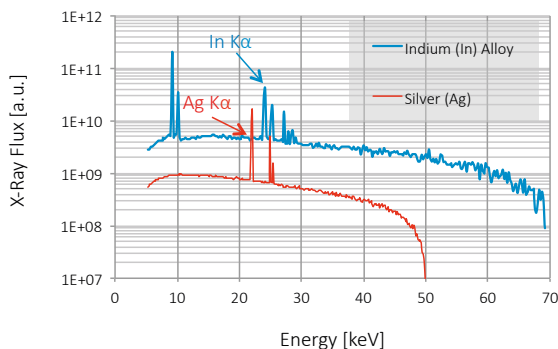


Spot Quality

Thanks to advanced electromagnetic focusing and correctional optics together with a high brightness LaB₆ cathode, a high quality e-beam focus is achieved. Together with a continuously generated smooth liquid target surface, the source produces X-ray spots of very high quality.

Line Profile Through Maximum

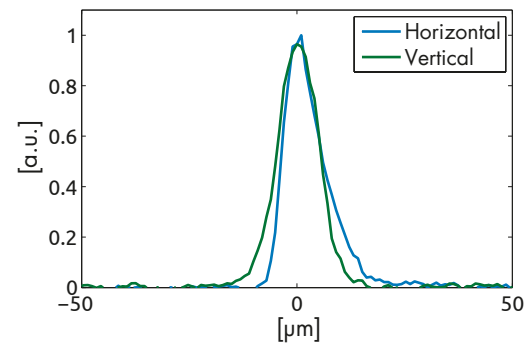

X-Ray Spot

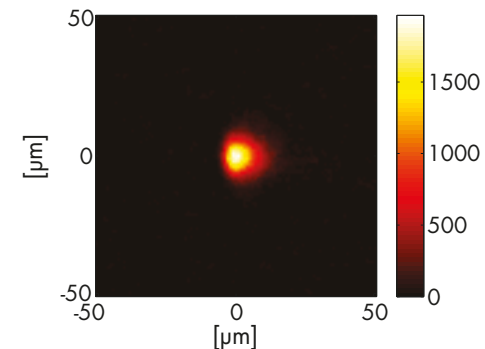


Tunable Size


Both the spot size and the aspect ratio can be tuned freely.

Source Stability

The spatial stability of the source is very high. The image below illustrates a spot centroid standard deviation of $< 0.1 \mu$ m over 24 hours, as taken with pinhole camera mechanically coupled to the source.

Spot Position Over 24 h

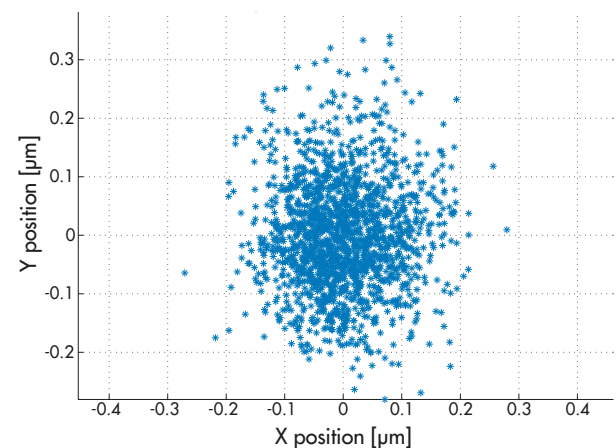



Excillum and its employees are dedicated to finding the best solution to our customers' needs. We work in close collaboration with our OEM and end customers to provide the optimum X-ray source for any application.

Please contact us with your specific requests!

Excillum AB
Jan Stenbecks Torg 17
164 40 Kista, Sweden

excillum

E-mail: sales@excillum.com
Website: www.excillum.com
Phone: +46-8-7509660

Although diligent care has been used to ensure that the information herein is correct, nothing contained herein should imply any representation or warranty as to the accuracy, currency or completeness of this information. The content hereof is subject to change without further notice. Please contact us for the latest version of this document or further information. Revision 2019-01-14.

Excillum, Metaljet, Metaljet D1, Metaljet D2, Metaljet D2+, Metaljet C2, NanoTube, and NanoTube N1 are registered trademarks or trademarks of Excillum AB. Excillum's X-ray sources are protected by several patents including, but not limited to, US Patents No. US 6 711 233, US 8 170 179, US 8 681 943, US 8 837 679, US 9 171 693, US 9 245 707, US 9 380 690, US 9 530 607, US 9 564 283, US 9 947 502, and Chinese Patents No. ZL 01816396.3, ZL 200780026317.0, ZL 200980155094.7, ZL 200980158566.4, ZL 201080070417.5, ZL 201180058633.2, ZL 201280075230.3, ZL 201410213235.9, ZL 201510020687.X, and other corresponding national patents and patent applications pending.